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Users are required to follow the safety and security advice.

Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

The cross roller bearing guides

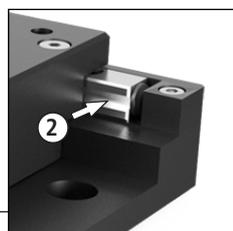
are lubricated by us with the **OWIS® SST.F2** grease.

Lubricating grease SST.F2 in a 5 ml applicator, part no. 90.999.0001

Relubrication should take place after about 500 operation hours, at least every 3 months.

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Lubrication procedure

1. Turn off the control unit and make sure that it cannot be turned on accidentally. It is then possible to move the carriage by hand.
2. Remove the used grease or dust particles from the guides with a dry, lint-free cloth. Cleansers such as benzine or petroleum are normally not necessary.
3. Insert the applicator cannula (1) into the open bore and fill in about 0.1 ml grease SST.F2.
4. Also apply a thin film of lubricating grease SST.F2 onto the raceways (2) of both crossed roller bearing guides.