## Goniometer Goniometers

## MOGO/TPM 150

/ 23.11.2020

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**Users are required to follow the safety and security advice.**

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Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

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The **worm gear** and the **guides** are lubricated by us with the **OWIS® SST.F8** grease.

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Lubricating grease SST.F8.L in a 5ml applicator, long, part no. 90.999.1008

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Relubrication should take place after about 700 operation hours, at least every 3 - 6 months. It is recommended to send the product to OWIS® for a complete service after 5 years.

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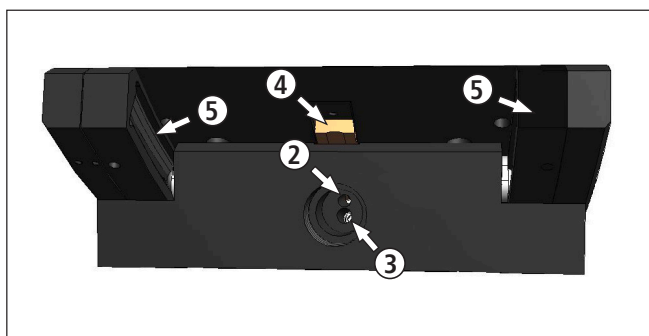
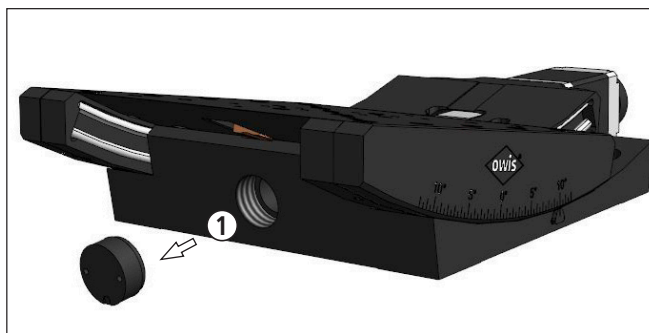
1. Move the MOGO slowly onto the limit switch (MAXDEC) and release it. Afterwards, turn off the control unit and make sure that it cannot be turned on accidentally.

2. Remove the screw (1) of the lateral bore, to uncover the maintenance bore (2) and the hexagon socket (3).

**Attention!** During the actual lubrication procedure the MOGO must be halted, so that the applicator needle cannot be captured and get screwed in by the worm.

3. Insert the applicator needle into the open maintenance bore up to the worm wheel segment (4) and fill in some grease SST.F8. Also apply a thin layer of SST.F8 evenly on the guides (5).

4. Move the goniometer manually by a quarter of its travel with a hexagon key (AF 3) and continue the process over the whole travel range.

5. Fix the screw (1) of the lateral bore.

6. To have the grease distributed equally, the axis should be moved over the whole travel two to three times.